

Mastering micropipetting

Micropipetting target practice

You will need (per pair of learners):

- P20 micropipette
- Tube of food dye
- Box of micropipette tips
- Laminated micropipetting target practice sheet
- Waste disposal container
- Non-permanent marker pen

1. In pairs, using a non-permanent marker pen, write your names on the lines on the laminated micropipetting target practice sheet.
2. Take it in turns to use the micropipette to accurately dispense the volumes of food dye shown into the centre of the circles on the laminated micropipetting target practice sheet.
3. Dispose of the used pipette tip in the designated waste disposal container.
4. Compare your results with your partner's to see if you have both mastered the skill of micropipetting!

Micropipetting target practice					
Name	20 µl	15 µl	10 µl	5 µl	2 µl
	○	○	○	○	○
	○	○	○	○	○